



Exceeding Your
Expectations
We Build
Trust

深圳市罗杰科技有限公司
SHENZHEN SCIFLOW TECHNOLOGIES CO LTD
www.sciflow.com eleflow@163.com

UD Cohesive Sampler



Unit Dose Sampler for Cohesive Powders

A popular sampler in the pharmaceutical industry for taking low volume (unit dose) cohesive powder samples. Allows accurate samples to be taken.

Type: Single Point Unit Dose

Material: 316 stainless steel

- Operation:
1. Insert the sampler into the product.
 2. At the required depth push down the handle and rotate it through 180°.
 3. Pull up sampler several millimetres (to fill the cell), then rotate handle 180° (to lock cell in closed position).
 4. Remove sampler

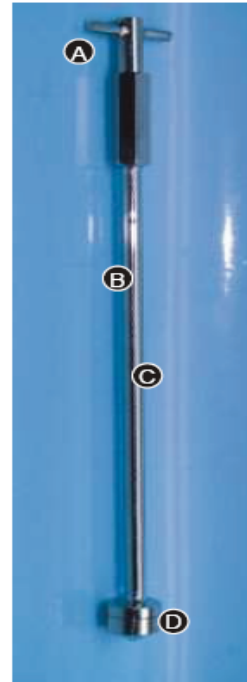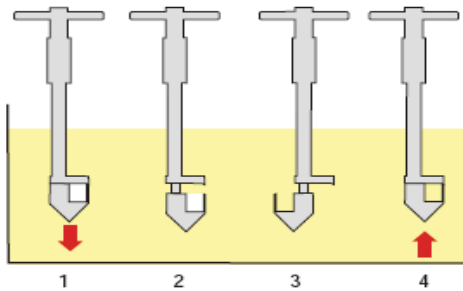


- (A) 'T' Bar Handle
- (B) Special lengths available on request
- (C) Polished finish
- (D) Sample Collecting Cell - different sized cells can be easily interchanged



'T'Bar Handle



Tip in open position

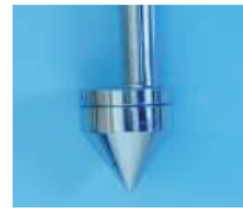

Tip in closed position



Part No.	Description
2000S-01	U-D Cohesive Sampler Set - consisting of: 1 off Handle, Sample Collecting Cells - 1 each of: 0.25ml, 0.5ml, 0.75ml and 1.0ml, 1 off Superior Sampler Case
2000A-1000	Sampler body and handle - 1000mm length
2000T-025	Sample Collecting Cell - 0.25ml
2000T-050	Sample Collecting Cell - 0.5ml
2000T-075	Sample Collecting Cell - 0.75ml
2000T-100	Sample Collecting Cell - 1.0ml
8001N-1100	Cleaning Brush - 1100mm length - narrow style
	Storage Cases - see page 79

深圳市南山高新区科苑南路留学生创业大厦 1406 Tel: 86-755-86329015/86329083 Fax 86-755-86329019
R1406 Overseas Chinese Scholar Venture Park, South Keyuan Road, Nanshan Dist., ShenZhen518057, China